

ANALAB®

P'Tips Cleaner

Pipette Tip Cleaner

*Better Sample Preparation
= Better Results!*



Convenient, Safe, and Effective Cleaning for Pipette Tips

P'Tips Cleaner removes contamination from pipette tips, readying them for trace metal determinations. Tips can be safely and simply cleaned with acid solutions using these gravity-fed devices.

- Contamination is leached from inside and outside of pipette tips
- Convenient batch cleaning of pipette tips
- Safe, closed system eliminates user contact with strong acids
- Cleaning performed on up to a full rack of tips at a time
- P'Tips cleaning process even allows tips to be reused if desired



**See it in
Action!**



How It Works

- Place pipette tips mounted in Cleaning Holder into P'Tips cleaning reservoir and replace lid
- Place PFA bottle containing cleaning solution in Fill Position – solution gravity feeds into reservoir and cleaning process begins
- After a period of time, simply move PFA bottle to Drain Position and solution drains out of reservoir
- Cleaning solution can be reused for several cycles and replaced as needed



Fill Position



Drain Position

Models

Two P'Tips Cleaner models are available to accommodate various pipette tip sizes.

P'Tips Model	Pipette Tip Sizes (μL)	Part Number
Medium	100 / 200 / 1,000	PT-M
XLarge	100 / 200 / 1,000 / 5,000 / 10,000	PT-XL



P'Tips Cleaner Medium


P'Tips Cleaner XLarge

Accessories

The P'Tips Holders and Transfer Racks are designed to match the configuration of commercial shipment racks that smaller pipette tips (200 and 1,000 μL) are commonly supplied in, which eliminates manual handling of the tips. Complete your P'Tips system by choosing the appropriate Holders and Transfer Racks to meet your laboratory's needs. High-purity precleaned pipette tips are also available.


P'Tips Holders

Supports tips during cleaning

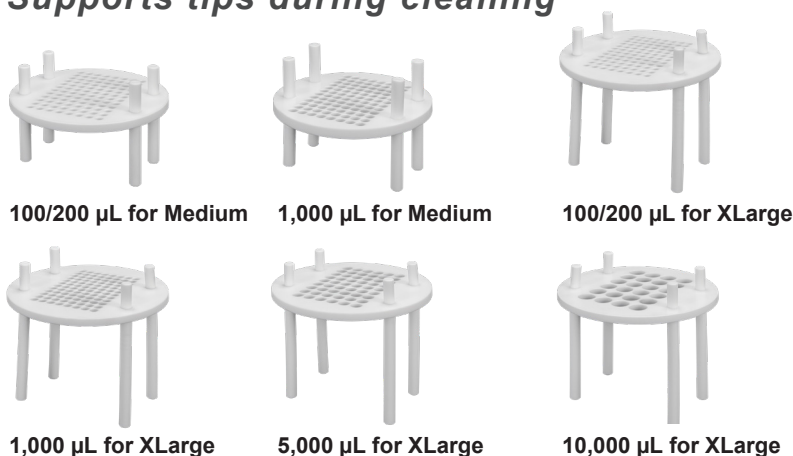

P'Tips Transfer Racks

Relocates tips in/out of Holder

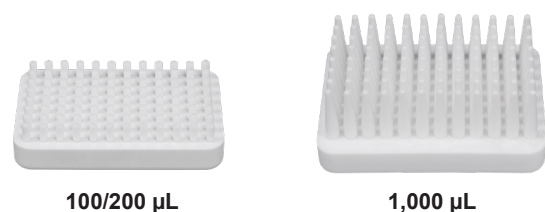

PFA Pipette Tips

High-purity precleaned



P'Tips Accessories	Part Number
100 / 200 μL P'Tips Holder for P'Tips Medium	PH-200-M
100 / 200 μL P'Tips Holder for P'Tips XLarge	PH-200-XL
1,000 μL P'Tips Holder for P'Tips Medium	PH-1000-M
1,000 μL P'Tips Holder for P'Tips XLarge	PH-1000-XL
5,000 μL P'Tips Holder for P'Tips XLarge	PH-5000-XL
10,000 μL P'Tips Holder for P'Tips XLarge	PH-10000-XL
100 / 200 μL P'Tips Transfer Rack	PTR-96-200
1,000 μL P'Tips Transfer Rack	PTR-96-1000

PFA Precleaned Pipette Tips	Qty.	Part Number
100 μL PFA Pipette Tips	10	ES-7000-0101
100 μL PFA Pipette Tips	96	ES-7000-0111
1,000 μL PFA Pipette Tips	10	ES-7000-1001
1,000 μL PFA Pipette Tips	60	ES-7000-1011